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- Ready to use biocidal product with broad spectrum
- Synergistic combination of 2 antimicrobial active ingredients
- Formaldehyde free
- Activity in compliance with European standards
- Made with WFI
- Double-bagged packaged
- Gamma irradiated

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### INDICATIONS

Cleaning and disinfection of floors, walls, surfaces and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

### COMPOSITION

Glutaraldehyde [CAS N°. 111-30-8 : 0,5 mg/g], didecylidiméthylammonium chloride [CAS N°. 7173-51-5: 1,2 mg/g].

### MICROBIOLOGICAL PROPERTIES

Bactericide in 5 min, 20°C (EN 1276, EN 13697).  
Yeasticide in 15 min, 20°C (EN 1650, EN 13697).  
Active in 5 min, 20°C against BVDV (surrogate of HCV) and PRV (surrogate of HBV).

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...): see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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## PHYSICO-CHEMICAL DATA

- Liquid colourless solution ready to use
- Soluble in all proportions in water
- Density at 20°C : approx. 1
- pH of product at 20°C : 3 to 5
- Non-corrosive towards metals (non oxidising).

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## INSTRUCTIONS FOR USE

To be used undiluted.  
Remove the first package before entering into the clean room. Remove the second package on first use.  
Apply a sufficient amount (+/- 30 ml/m<sup>2</sup>) respecting the indicated contact time.  
If necessary, remove the excess after the contact time.  
For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

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## PRECAUTIONS FOR USE

Dangerous - Please follow the instructions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products).  
Use the biocides with precaution. Before use, read the label and the product information.  
Store between +5°C and +35°C.  
Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

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